


MELING
BIOLOGY & MEDICAL

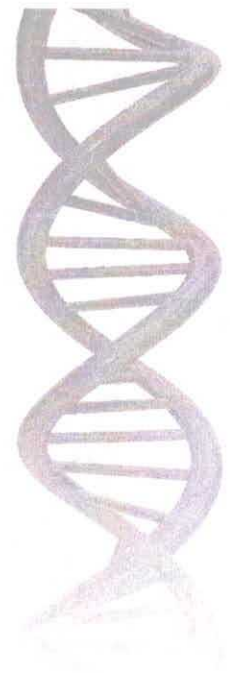

Operation Manual

The following product models are applicable:

Biomedical Freezer

DW-YL450

Ultra Low Temperature Freezer

DW-FL450

Combined Refrigerator and Freezer

YCD-EL450 / YCD-FL450

Zhongke Meiling Cryogenics Company Limited

MELING
BIOLOGY & MEDICAL












Better science, Better life!










www.melingbiomedical.com

zkmeiling@zkmeiling.com

Table of Contents

1. Application Notes	1
2. Safety Instructions	2
3. Precautions in Use	4
4. Product Installation	4
4.1 Installation Environment	4
4.2 Installation Site	4
4.3 Preparation before Use	5
4.4 First Power-on	5
4.5 Operation after Power Failure	5
5. Product Composition and Overview	6
6. Operating Instructions	7
6.1 Function Introduction	7
6.2 Optional Function	8
7. Defrost , Discontinuation and Maintenance	10
7.1 Equipment Maintenance	10
7.2 Equipment Discontinuation	10
7.3 Maintenance, Replacement, and Recovery of Rechargeable Batteries	10
8. Troubleshooting and Maintenance Services	11
9. Specifications	12
10. Packing List	12
11. Electrical Schematic Diagram	13

	Do not place the equipment in areas exposed to the sun or rain, so as to prevent danger such as short circuit or overheating.
	Do not tilt or lay the equipment sideways, and do not impact the equipment body; Refrigeration systems are installed in the equipment, which is easy to be damaged by tilt or impact.
	Please place the equipment in a dry and dust-free environment to avoid risks such as overheating, and short circuit.
	In case of unexpected sound, smell, smoke, etc. when the power is turned on, please unplug the power in time and contact the manufacturer or supplier.
	Please place the equipment in a dry and ventilated environment, and ensure that the equipment vents and instrument surfaces are not blocked or shielded by walls or other objects; Do not use it in a poorly ventilated environment to prevent damage caused by heat released by equipment.
	It is forbidden to disassemble and modify this equipment without authorization, so as to avoid potential safety hazards. In this case, MELING BIOLOGY & MEDICAL will not bear any responsibility for quality accidents.
	It is forbidden to put inflammable and explosive dangerous goods, strong corrosive acids, alkalis and other items unsuitable for the equipment in the equipment.
	When storing toxic, harmful or radioactive materials, please use the equipment in safe areas. Improper use may cause harm to human health or environment.
	Metal objects such as nails or iron wires shall not be inserted into any aperture and gap or any outlet of the equipment, otherwise electric shock or injury may be caused due to accidental contact between the above objects and moving parts.
	In order to ensure the normal operation, ventilation and heat dissipation of the equipment, the back, left and right sides of the freezer body should be at least 30 cm away from the wall, and the air inlet and air outlet shall not be blocked by obstacles!
	This equipment must be connected to a ground wire.

	Note: Failure to observe the precautions may result in personal injury or equipment failure and related property losses.
	It is forbidden to store living animals, flowers or other items with strict temperature requirements in the equipment.
	When the equipment is running, do not touch the inner surface of the cabinet without wearing protective gear.
	Hold the handle and close the door to avoid pinching your fingers; When the equipment is not used for a long time, please unplug it and pack it for storage.
	When restarting the equipment after power failure or power off, please check the equipment settings first, otherwise the stored items may be damaged due to the change of settings.
	The equipment can be used for item preservation, not as production equipment!
	Keep the keys properly, so as to avoid accidents when children open the door accidentally.
	When handling the equipment, please be careful not to tip over the equipment, so as to prevent equipment damage or personal injury.
	When handling, it shall be lifted from the bottom, with the inclined plane not be greater than 45°, and it shall be handled with care. Please use the equipment in safe areas. Improper use may cause harm to human health or environment.

4.3 Preparation before Use:

1. Remove the outer packaging of all products (including the protective foam in the packaging box)



Warning: Don't put the plastic bags within the reach of children, so as to prevent suffocation accidents.

2. Inventory of accessories: Please check the accessories and materials according to the packing list.
3. Cleaning: Clean the product once before use.

4.4 First Power-on

When using the equipment for the first time, please follow these steps:

1. After the equipment is placed, leveled and cleaned, it shall stand for more than 24 hours, and then power it on to ensure the normal operation of the equipment.
2. Under no-load condition, connect the power cord to a special outlet with appropriate specifications.
3. After powering on, turn on the power switch of the equipment.
4. Check whether the operating temperature of the equipment reaches the required value, observe the normal start and stop of the equipment for more than 24 hours, and put a small amount of items in the freezer after confirming the normal performance.
5. Please store items in batches, with the items not exceeding 1/3 of the freezer volume each time. Ensure that the equipment is running properly after shutdown for more than 12 hours before putting in the next batch of items).
6. Try not to open the door during cooling, otherwise the temperature will rise.

4.5 Operation after Power Failure

1. The equipment has a memory function for the set value. When the power is restored after power failure, the equipment will continue to operate before power failure.
2. Once the equipment is powered off, it requires 5 minutes before it can be switched on again, so as to avoid damaging the compressor.
3. We guarantee the normal operation of this equipment under certain conditions, but we are not responsible for any loss or damage of stored items after power failure.



Notes:

- ◆ The transportation and movement of equipment is realized by casters.
- ◆ A special person shall be responsible for checking and recording the running status of the equipment every day (record and check once every 2-4 hours). In case of failure or shutdown, the temperature in the freezer will rise. If it cannot be repaired in a short time, please take out the stored items and transfer them to a place that meets the temperature requirements for storage to avoid damage to the items.
- ◆ Before putting items into the equipment, it shall be confirmed in advance whether the temperature range of the equipment meets the temperature requirement of the items, so as to avoid damage to stored items due to the difference between the settable temperature of the equipment and the required temperature of the items. Please pay attention not to block the air outlet and air inlet when putting items into the equipment.
- ◆ Due to the refrigeration inertia, there is a certain difference between the actual display temperature and the set temperature of the equipment, which is a normal phenomenon.
- ◆ The equipment is an item storage equipment, which cannot be used for routine production operations. It is strictly forbidden to put too many items which are relatively hot into the equipment at one time, otherwise the compressor will run for a long time, and be burned due to high temperature. Items must be put in batches, so as to ensure that the freezer is cooled step by step until the temperature required for storing items is reached.
- ◆ Electrical appliances without production license shall not be used inside the equipment.
- ◆ Do not change the set temperature frequently in a short time, otherwise the expected setting effect may not be achieved due to the large temperature inertia; Ensure that there is a certain air circulation space around the cabinet when putting in items, especially do not block the temperature sensor in the cabinet (for collecting the temperature of the cabinet), otherwise it will affect the stability and accurate control of the temperature in the freezer body!
- ◆ Items shall not be placed directly at the bottom of the freezer, but on the bottom shelf, otherwise the refrigeration effect of the equipment will be affected.
- ◆ When putting in items, if the moisture content of the items is too much or too little, it will affect the humidity change in the cabinet, so it is best to keep the items sealed; The humidity of the working environment will affect the change of humidity in the cabinet, especially if the door is opened too frequently and the door is not closed properly.



Warning: Children are not allowed to play with this equipment as a game prop, otherwise the injury or loss caused therefrom will be at their own risk.

6.2 Function Settings

1. After powering on, the machine will enter into the working state;

2. User parameter settings:

Unlock: Under normal operation, press **▲** and **▼** key for 3 seconds at the same time, and the digital tube in the upper chamber will display "PS1" and the digital tube in the lower chamber will display the parameter code "0000." Press **▲**, enter the password "0005" (when entering the user menu password, enter "0099" to restore the key lock password to the default "0005."), and unlock it. Unlock and press **■** key for 3s, and the digital tube will display the parameter code "PS1," and enter the setting and adjustment parameters.

① Use **■** to scroll the parameters, and the display order is as follows: b1→b2→Set1→H01→L01→Set2→H02→L02→n→y→r→S→F→Pt→tH1→P1→Ps1.

② Use **▲** or **▼** to increase or decrease the value;

③ Use **■** to temporarily store the modified values and return to the display parameters;

④ If other parameters are modified, repeat steps ①~③.

⑤ Press **■** for more than 3s, save the modified parameters and return to the display parameter category.

3. Press **■** for more than 3s, or press no key in 60S to exit the parameter setting program.

4. Parameter display

No.	Menu item	Parameter Range	Suggested settings	Remarks
1	b1	---	---	Hardware version
2	b2	---	---	Software version
3	Set1	DW-YL series: -10~-25 DW-FL series: -20~-40 YCD-EL series: -10~-26 YCD-FL series: -10~-40	DW-YL series: -25 DW-FL series: -40 YCD-EL series: 5 YCD-FL series: 5	Upper chamber temperature setting
4	H01	0. 0~10. 0	5. 0	Set value of high temperature alarm set+H; When H =0, High temp alarm is disabled; When the alarm is over high temp alarm set, H1 will be displayed on the controller
5	L01	0. 0~10. 0	5. 0	Set value of low temperature alarm set-L; When L =0, Low temp alarm is disabled; When the alarm is below low temp alarm set, L1 will be displayed on the controller
6	Set2	DW-YL series: -10~-25 DW-FL series: -20~-40 YCD-EL/FL series: 2~8	DW-YL series: -25 DW-FL series: -40 YCD-EL series: -20 YCD-FL series: -40	Lower chamber temperature setting
7	H02	0. 0~10. 0	5. 0	Set value of high temperature alarm set+H; When H=0, cancel the alarm; When the temperature is too high, the high temperature alarm displays H02
8	L02	0. 0~10. 0	5. 0	Set value of low temperature alarm set-L; When L=0, cancel the alarm; When the temperature is too high, the high temperature alarm displays L02
9	Pt	0-240 min	20	Print interval
10	tH1	20.0~50. 0℃	50	Upper limit of ambient temperature alarm
11	P1	Automatic heating mode 1 Automatic heating mode 23. Automatic heating mode 34. Always on5. Always off	1.(Set to 4 when the door is with condensation)	Mode 1: It shall be judged as once after the door is opened and closed once and heated for 5min (time setting). If the door is opened and closed again during the heating period, the heating time will be updated again; Mode 2: After the compressor is running. Heater on; When the compressor stops, the heater is delayed for 1 minute for turn-off; Mode 3: When the humidity in the cabinet is more than 80%, the door heating is on, and when the humidity in the cabinet is moderately less than 60%, the door heating is off; Mode 4: Door heating is always on; Mode 5: Door heating is always off.
12	Ps1	0000~9999	0005	User menu password settings

5. Alarm display

Alarm Code	Default Description
H01	Upper chamber high temperature alarm
L01	Upper chamber low temperature alarm
H02	Lower chamber high temperature alarm
L02	Lower chamber low temperature alarm
H03	High ambient temperature alarm
door	Door opening alarm
PF	Power failure alarm
bL	Low battery alarm
EE	Communication failure
E2	Upper chamber temperature sensor alarm
E4	Lower chamber temperature sensor alarm
Er	Recorder not connected
Pr	USB Data logging failure
LoF	Recorder not started

6.3 Optional Function

1. Printer

Automatic printing:

When the printing interval P_t is not 0, the printer prints the current data every P_t minutes.

Note: When the printing interval P_t is less than the recording interval SC_y , it is printed according to the recording interval; the remaining printing intervals are printed as an integral multiple of the recording interval SC_y , that is: the printing interval P_t should be set as an integral multiple of the recording interval.


Manual printing:

When the key is unlocked and the working voltage is normal, press the print key to display "P01" by the digital tube, press the up key or down key to adjust "P00~P07." Press the set/mute multiplexing key or print key, and the printer will cancel the manual printing (P00) or print the data recorded in the previous days (1 ~ 7), with the data interval same as the automatic printing interval.

Printing paper installation:

The printing paper has been installed at the factory. When the paper is used up after a long time, you may purchase the same paper for installation. The installation steps are as follows:

- ① Click the cylindrical button on the printer to open the cover of the printer;
- ② Put the printing paper into the printer box and pull the paper head of the roll so that the paper slightly exposes from the upper port of the box cover after the box cover of the printer is closed;
- ③ Close the box cover.

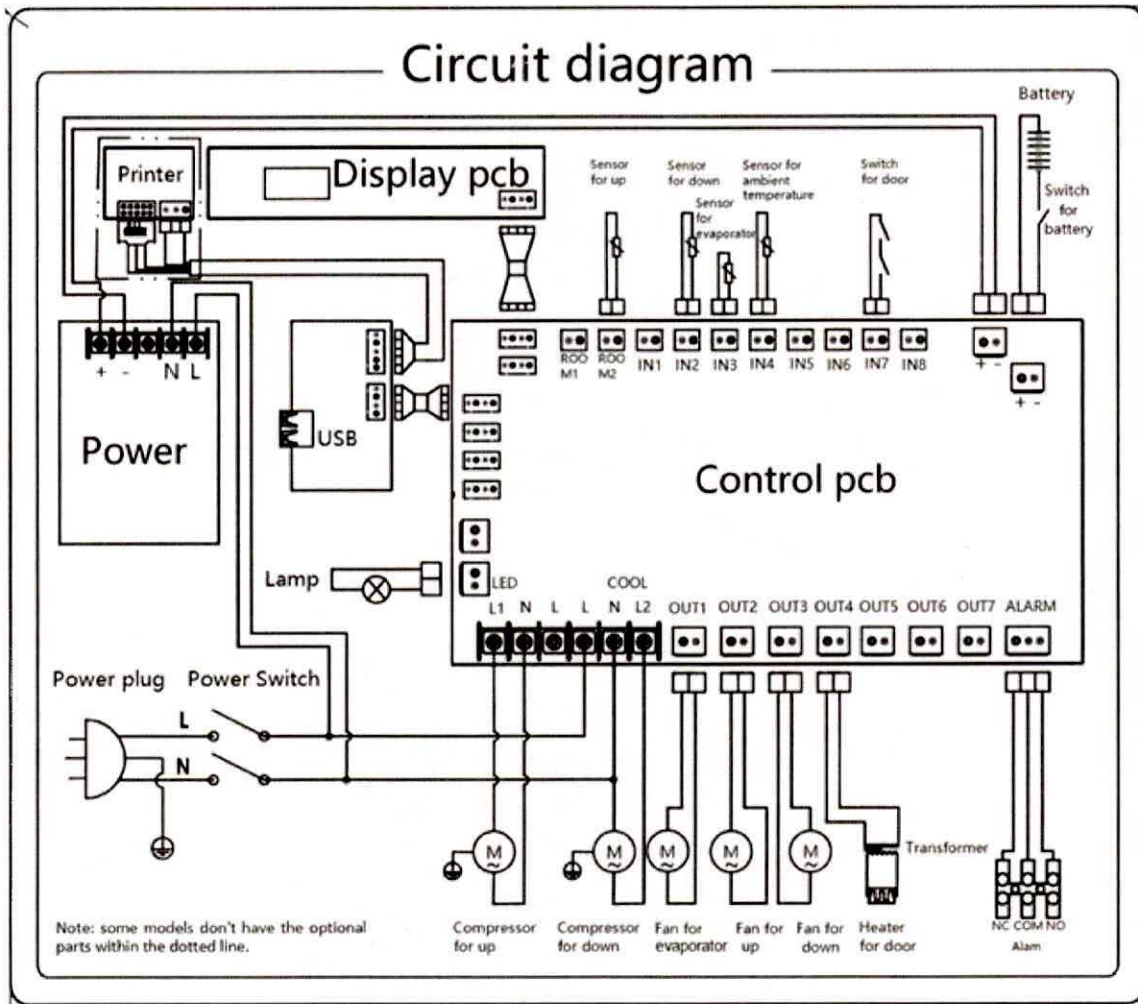

YCD-EL450/YCD-FL450

* The technical data and electrical schematic diagram shall be subject to the nameplate on the product and the electrical schematic diagram in case of any improvement of the product design. We sincerely apologize for any inconvenience that might be caused.



Zhongke Meiling Cryogenics Company Limited
Address: No 1862 Zishi Road, Hefei City, Anhui, P.R. China
Production Address: No.1862 Zishi Road, Economic and
Technological Development Zone, Hefei City
Post Code: 230601
Email: zkmeiling@zkmeiling.com; technical.service@zkmeiling.com
Website: www.melingbiomedical.com
Material Code:890305660
Production Date: See nameplate on the freezer body
Prepared in: June 2020